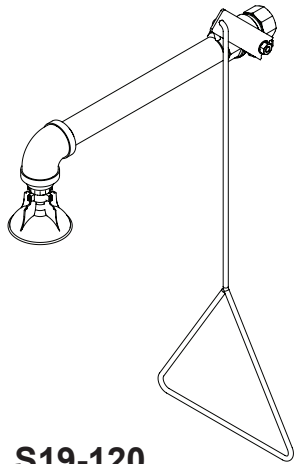
# Installation

## S19-120

Drench Shower Horizontal Supply  
with Plastic Showerhead

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
S19-120

## S19-120A

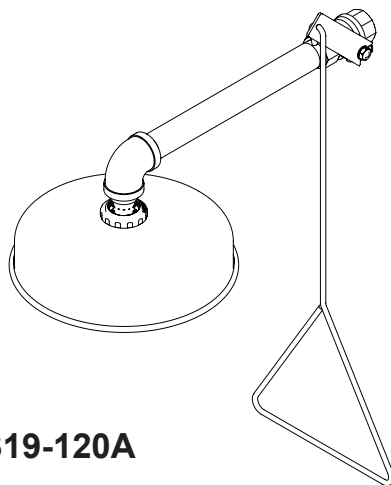Drench Shower Horizontal Supply  
with Plastic Showerhead and  
Stainless Steel Shroud

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S19-120A

Read the instructions in this manual before beginning installation. Save these instructions and refer to them for inspection, maintenance and troubleshooting information.

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## Table of Contents

|   |   |
|---|---|
| Safety Information .....                              | 3 |
| Supplies Required.....                                | 4 |
| Assemble components.....                              | 4 |
| Connect water supply.....                             | 4 |
| Assembly of Components and Parts List - S19-120.....  | 5 |
| Assembly of Components and Parts List - S19-120A..... | 6 |

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For questions regarding the operation, installation or maintenance of this product, visit [bradleycorp.com](http://bradleycorp.com) or call 800.BRADLEY (800.272.3539).

Product warranties and parts information may also be found under "Resources" on our website at [bradleycorp.com](http://bradleycorp.com).

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### To ensure proper operation:

#### Installation

This fixture has been designed and 3rd party tested to meet ANSI/ISEA Z358.1. The system must be tested periodically to be sure it is functioning properly to meet the standard. Please refer to the ANSI/ISEA Z358.1 standard on testing requirements.

The installation, maintenance, and location of all safety drench showers, eye, and eye/face washes must comply with the requirements of ANSI/ISEA Z358.1. Workers who may come in contact with potentially hazardous materials should be trained regarding the location and proper operation of emergency equipment per ANSI/ISEA Z358.1.

ANSI/ISEA Z358.1 requires this unit to be used with a clean, potable, uninterrupted supply of tepid water. Bradley plumbed emergency fixtures require a minimum of 30 psi flowing pressure in order to function as designed and meet the ANSI/ISEA Z358.1 standard.

Before installation make sure that this equipment will meet the requirements of the potential hazardous contaminants in your location.

Installation of this system must be completed by a qualified plumber in compliance with all national and local codes. Compliance and conformity to drain requirements and other local codes and ordinances is the responsibility of the installer.

#### Weekly Inspections

To ensure all emergency shower and eye/face wash fixtures are ready to be used when needed, regular checks and maintenance must be conducted. Weekly inspections must be conducted on all safety equipment to make sure a suitable flushing fluid supply is present and to make sure the equipment is in good operating condition.

#### Water Temperature

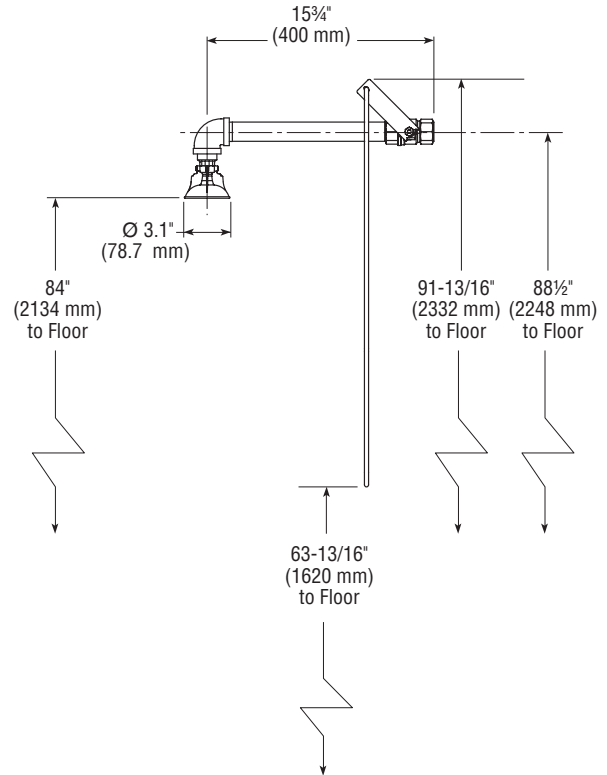
ANSI/ISEA Z358.1 requires tepid water. Suitable range is 60°F to 100°F (16°C to 38°C). Personal injury is possible outside this temperature range.

## Supplies Required

- Teflon tape and pipe sealant
- Piping to 1" NPT water supply inlet
- Adequate supply pipe supports
- Minimum 4" (102 mm) drain to accommodate 30 gallons (115 liters) per minute discharge for shower waste
- Sign-mounting hardware

## Assemble components

1. Assemble the components as shown on page 4 (S19-120) or page 5 (S19-120A).
2. Apply pipe sealant to all male-threaded pipe joints.
3. Use the rubber grip pad provided or a strap wrench around pipes when tightening to prevent marring. Place the grip pad on the pipe, then put the wrench over the grip pad and turn the pipe with the wrench.
4. Bottom edge of showerhead should be 84" (2134 mm) from the floor.

S19-120


## Connect water supply

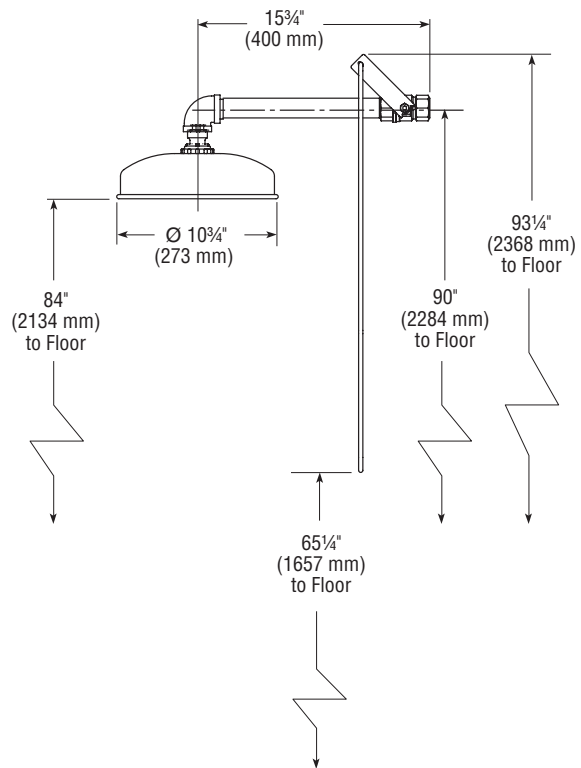**NOTICE! Do not rely on Bradley's Drench Shower to support supply piping.**

1. Connect water supply piping to 1" NPT inlet on unit (piping supplied by installer). Provide adequate supports (supplied by installer) for supply pipe using pipe hangers or other means.
2. Mount the safety sign to the wall using sign-mounting hardware supplied by installer.
3. Open water supply lines. Test for leaks and adequate water flow.


All dimensions assume standard thread engagement. Variations in manufacturing allow for +/- 1/8" (3 mm) per threaded joint. To find the tolerance of a dimension, add the number of thread joints in between a dimension and multiply it by 1/8" (3 mm).

S19-120A

# Assembly of Components and Parts List - S19-120

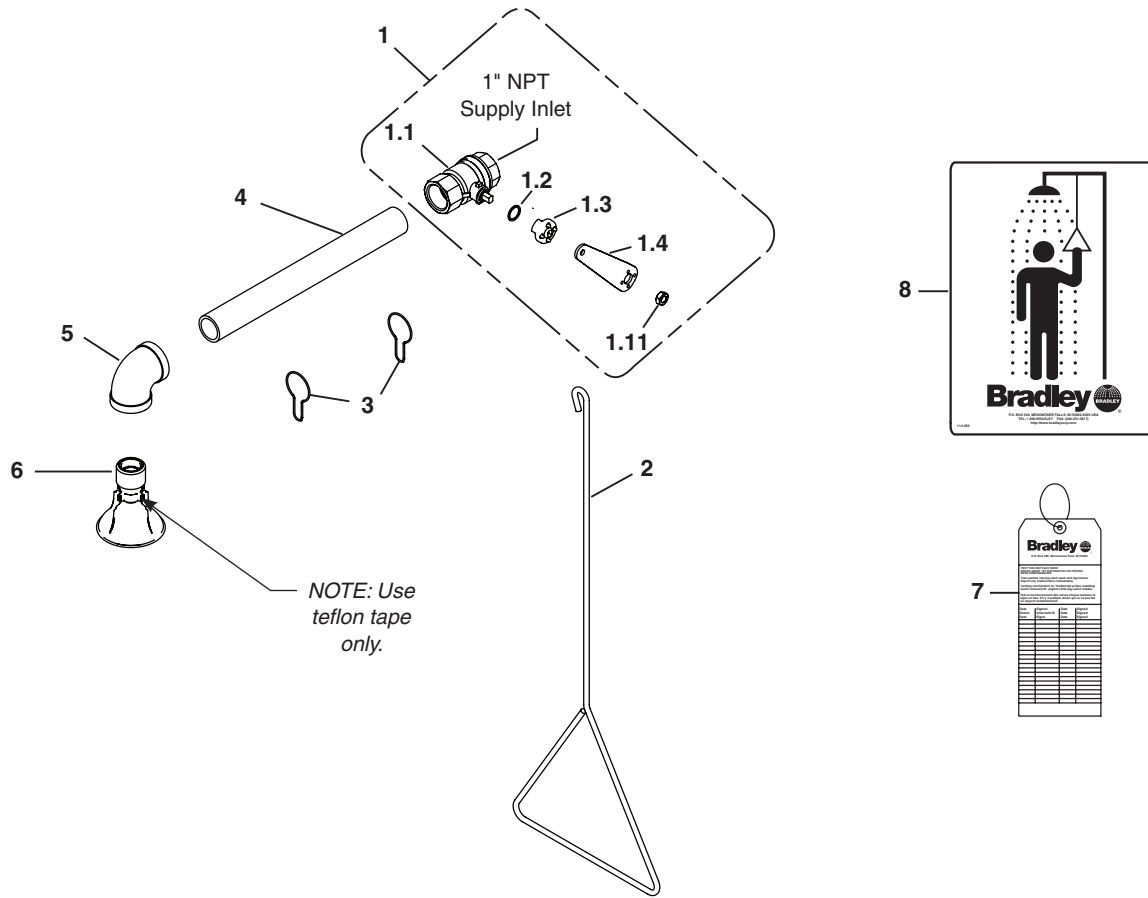
NOTE: Items 1.1–1.5 come preassembled as Item 1.

| Item | Part No. | Qty. | Description                   |
|------|----------|------|-------------------------------|
| 1    | S30-059  | 1    | Stay-Open Ball Valve Assembly |
| 1.1  | S27-278  | 1    | Ball Valve with Jam Nut       |
| 1.11 | 110-214  | 1    | Jam Nut (only)                |
| 1.2  | 124-048  | 1    | Washer                        |
| 1.3  | 153-489  | 1    | Adapter                       |
| 1.4  | 128-201  | 1    | Handle                        |
| 2    | 128-156A | 1    | Pull Rod                      |

| Item | Part No.  | Qty. | Description                 |
|------|-----------|------|-----------------------------|
| 3    | 151-001   | 2    | Curtain Hook                |
| 4    | 113-006LK | 1    | Pipe 1" NPT x 12" Long      |
| 5    | 169-847   | 1    | 90° Elbow 1" NPT            |
| 6    | S24-188   | 1    | Plastic Showerhead Assembly |
| 7    | 204-421   | 1    | Emergency Tag               |
| 8    | 114-050   | 1    | Safety Sign                 |
| 9    | 269-915   | 1    | Grip Pad (not shown)        |

# Assembly of Components and Parts List - S19-120A

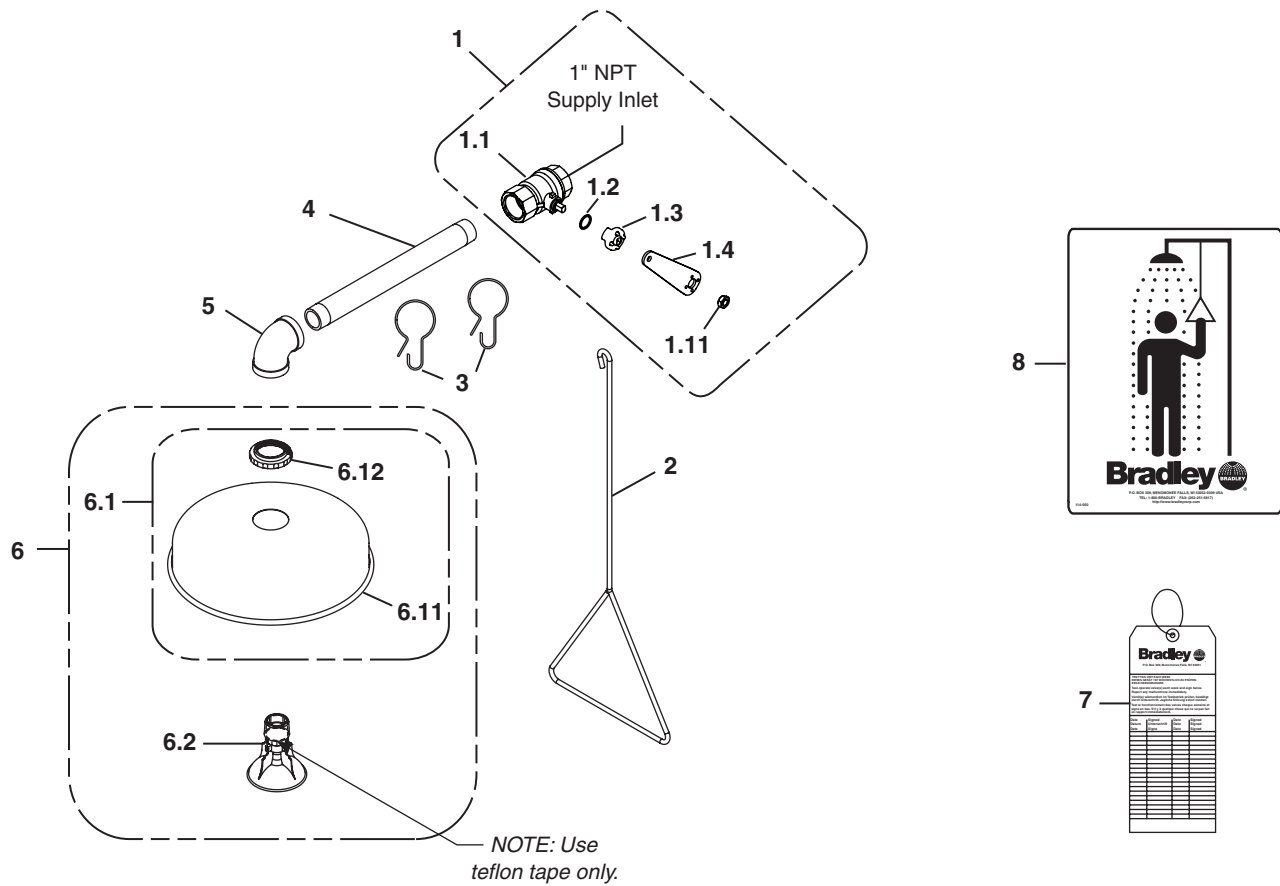NOTE: Items 1.1–1.5 come preassembled as Item 1.  
Items 6.1–6.2 come preassembled as Item 6.

| Item | Part No.  | Qty. | Description                   |
|------|-----------|------|-------------------------------|
| 1    | S30-059   | 1    | Stay-Open Ball Valve Assembly |
| 1.1  | S27-278   | 1    | Ball Valve with Jam Nut       |
| 1.11 | 110-214   | 1    | Jam Nut (only)                |
| 1.2  | 124-048   | 1    | Washer                        |
| 1.3  | 153-489   | 1    | Adapter                       |
| 1.4  | 128-201   | 1    | Handle                        |
| 2    | 128-156A  | 1    | Pull Rod                      |
| 3    | 151-001   | 2    | Curtain Hook                  |
| 4    | 113-006LK | 1    | Pipe 1" NPT x 12" Long        |
| 5    | 169-847   | 1    | 90° Elbow 1" NPT              |

| Item | Part No. | Qty. | Description                      |
|------|----------|------|----------------------------------|
| 6    | S24-189  | 1    | Stainless Steel Showerhead Assy. |
| 6.1  | S24-193  | 1    | Stainless Steel Shroud Kit       |
| 6.11 | 187-053  | 1    | Showerhead Shell                 |
| 6.12 | 154-147  | 1    | Ring                             |
| 6.2  | S24-188  | 1    | Plastic Showerhead Assembly      |
| 7    | 204-421  | 1    | Emergency Tag                    |
| 8    | 114-050  | 1    | Safety Sign                      |
| 9    | 269-915  | 1    | Grip Pad                         |

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